

# Loading dye (10X)

## Contents

15 %     Ficoll 400  
0.25 %   Bromophenol blue

### For 50 ml

7.5 g     Ficoll 400  
0.125 g   Bromophenol blue  
→ to 50 ml with dH<sub>2</sub>O  
→ store at 4°C

Ficoll 400 - Fisher BioReagents: BP525

Bromophenol blue - Acros Organics (via Fisher): AC15134-0250

### NOTES:

Make up in 50 ml centrifuge tube. Will be difficult to get into solution - leave O/N on orbital shaker.